



Alpha and Beta separation results with Triathler

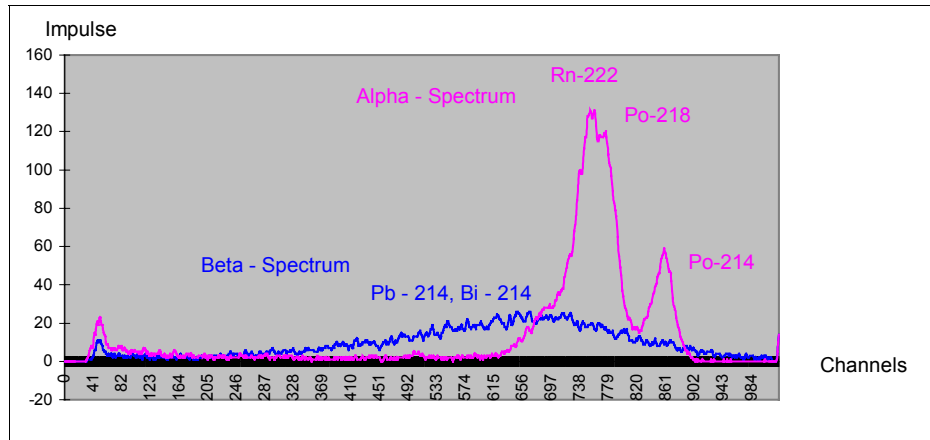

The 2-dimensional plot of alpha- and beta-spectra of Rn 222

Alpha – Beta separation is performed by the method of pulse length index; long alpha pulses are measured in the alpha channel, the significant shorter beta – pulses in the beta channel. The above figure (above) shows a typical Rn-222 spectrum, performed by Triathler™. In the 2-D spectrum (down) x – axis is energy, y–axis is pulse length and z – axis is intensity. The horizontal line is the pulse length index PLI, which discriminates alpha and beta pules, the vertical lines

correspond with the windows settings for alphas (top) and beta windows settings (down). PLI and windows can be set individually in order to perform best alpha - beta separation settings. The tool of the 2-D picture with Triathler™ Software is automatically created by the help of Excel –Macros. This tool enables optimal settings for Rn-222, U-234/U-234 or a - wipe test applications.

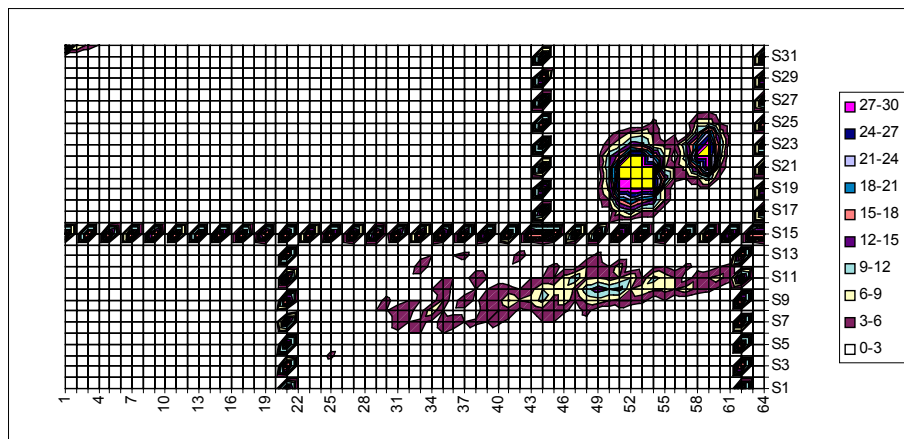

Product Information

Finland and International Hidex Oy tel +358-2-2750557 fax +358-2-2410075 info@hidex.com **Canada** Gamble Technologies Limited Tel: 905 812 9200 info@gtl.ca **France** ScienceTec Tel +3316446 2400 info@sciencetec.fr **Germany (Austria, Czech Republic)** FCI Frenzel Consulting & Instruments Tel +49 7082 948 290 eberhardfrenzel@aol.com **Hong Kong** Line Analytics Life Sciences Ltd. Tel + 852 2578 5839 line@lineanalytics.com **Japan** Sowa Trading Co Inc. Tel +81 3 3862 2700 tokyo@sowa-trading.co.jp **Korea** SIMS Co.Ltd Tel +82 2 576 9727 dmlee@sims.com **Russia** Pribori Oy c/o Wallac Oy Tel +358 2 267 8111 maria.kivisto@pribori.pp.fi **Scandinavia:** Reaction-Lab Oy Tel +358 2 410 1746 info@reaction-lab.com **Spain** Izasa s.a. Tel +34 91 6630500 jbruna@izasa.es **Switzerland** Uehlinger Pfiffner AG Tel +41 62 7212281 mail@uehlinger-pfiffner.ch **United Kingdom** Southern Scientific Ltd. Tel +44 1903 6044000 info@southern-scientific.co.uk **U.S.A** Bioscan Inc Tel + 1 202 338 0974 sales@bioscan.com

www.hidex.com